

## POLARIZATION-MAINTAINING FIBERS

### WHAT MAKES PM FIBERS MAINTAIN THE POLARIZATION?

In [polarization-maintaining single-mode fibers](#) (PM fibers), the fiber symmetry is broken by integrating stress elements in the fiber cladding. The light is then guided in two perpendicular principle states of polarization with different propagation constants – the fast and the slow axis. The linear polarization of light coupled into one of these axes is maintained. The linearly polarized laser radiation is conventionally coupled into the slow axis because of its lower sensitivity to fiber bending.


If light is coupled into both axis then the resulting polarization is elliptical (if the coherence length of the source is larger than the phase difference). Strain and temperature variations, however, change this arbitrary elliptical state. Thus it is important to exactly align the polarization axis of the laser source with the polarization axis of the fiber e.g. using the [Polarization Analyzer SK010PA](#).

Different types of polarization-maintaining fibers are designed depending on the geometry of the stress elements: "PANDA" fibers, "Bow-Tie" fibers or "Oval-Inner Clad" fibers.

The polarization-maintaining fiber cables made by Schäfter+ Kirchhoff typically use fibers of type PANDA. The slow axis is aligned with the index key of the FC type fiber connector with high precision ( $<1.5^\circ$ ).


The fiber cables made by Schäfter+Kirchhoff typically have a [polarization extinction](#)  $\geq 200:1$  (23 dB) or  $\geq 400:1$  (26 dB) for  $\lambda \geq 780$  nm (except for fibers end caps).

The fiber cables with [end caps](#) made by Schäfter+Kirchhoff typically have a polarization extinction  $\geq 200:1$  (21 dB) or  $\geq 400:1$  (23 dB) for  $\lambda \geq 780$  nm.

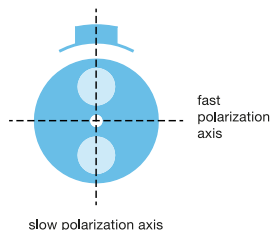


Polarization-maintaining single-mode fibers guide coupled radiation in two perpendicular principle states, the fiber polarization axes (also called the slow and fast axis, see figure on the left).

*Polarization axes of a PM fiber*

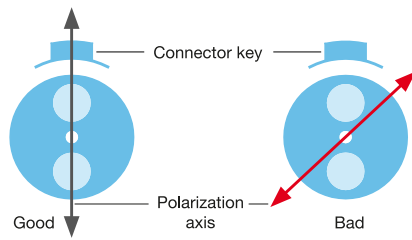

Axis orientation of a PM fiber

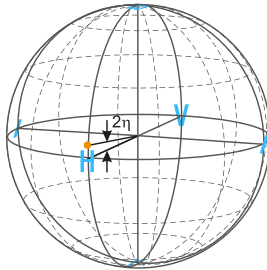

### COUPLING INTO THE FAST AXIS OF A PM FIBER

Conventionally the linearly polarized laser radiation is coupled into the slow axis because of its lower sensitivity to fiber bending.

You can also couple light into the fast axis of a PM fiber cable. However, the bandwidth of the fiber is severely reduced and it is only advisable to couple in a wavelength close to the cut-off.

Please also keep in mind that Schäfter+Kirchhoff only specifies and tests as well as supplies measurement data for radiation coupled into the slow fiber axis. Measurement and test result will vary for coupling into the fast axis.



SK010PA

The [polarization extinction ratio PER](#) of fiber-coupled radiation is the ratio between the optical power levels coupled to the two polarization axes of the fiber. The [Polarization Analyzer SK010PA](#) is used to optimize the polarization alignment of the polarization axis of the light source to the polarization axis of the fiber by rotating the input polarization axis of the source, see figure on the left.

- Good alignment: If the connector index key is in axis with the slow axis of the fiber and the laser polarization axis the alignment is good
- Bad alignment: If the connector index key is in axis with the slow axis of the fiber but the laser polarization axis is rotated the alignment is bad and the resulting output beam is arbitrarily elliptically polarized

### CHARACTERIZING POLARIZATION-MAINTAINING FIBERS (PM FIBERS)

More information on how Schäfter+Kirchhoff characterizes PM fibers can be found [here](#).

### POLARIZATION ANALYZER

More information of the application of the Polarization Analyzer can be found in the [Technotes](#) section as well as on the [product page](#). There is also an [article](#) that gives a good overview.

This is a printout of the page <https://sukhamburg.com/support/technotes/fiberoptics/cablebasics/pmc.html> from 6/15/2025

## CONTACT

For more information please contact:

Schäfter + Kirchhoff GmbH

Kieler Str. 212

22525 Hamburg

Germany

Tel: +49 40 85 39 97-0

Fax: +49 40 85 39 97-79

[info@sukhamburg.com](mailto:info@sukhamburg.com)

[www.sukhamburg.com](http://www.sukhamburg.com)

## LEGAL NOTICE

**Copyright 2020 Schäfter+Kirchhoff GmbH. All rights reserved.**

Text, image, graphic, sound, video and animation files and their arrangement on Schäfter+Kirchhoff GmbH webpages are protected by copyright and other protective laws. The content may not be copied for commercial use or reproduced, modified or used on other websites. [\[more\]](#)