
### FARADAY ISOLATOR: WORKING PRINCIPLE

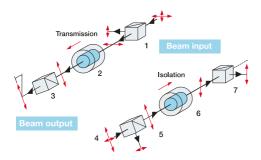

Faraday isolator scheme

## Effects on the input beam

- 1 Polarizing beam splitter eliminates the s-polarized portion (typically 1%) of the laser (diode) radiation
- 2 Faraday crystal in a strong magnetic field rotates the input polarization plane 45°
- 3 Polarizing beam splitter exactly aligned with the output polarization plane, i.e. by 45°with respect to the input polarization

### Effects on the back-reflected beam

- 4 Reflective or scattering surfaces cause polarized or depolarized reflected signals
- 5 Polarizing beam splitter (3) eliminates the s-polarized portion of the reflected signal
- 6 Faraday rotator (2) rotates the polarization plane by a further 45° so that the polarization plane of the reflected signal is now rotated by 90° from the polarization plane of the input beam
- 7 Polarizing beam splitter (1) blocks the reflected signal

This is a printout of the page <a href="https://sukhamburg.com/support/technotes/fiberoptics/multicube/48fi.html">https://sukhamburg.com/support/technotes/fiberoptics/multicube/48fi.html</a> from 7/6/2025

## CONTACT

For more information please contact: Schäfter + Kirchhoff GmbH Kieler Str. 212 22525 Hamburg Germany

Tel: +49 40 85 39 97-0 Fax: +49 40 85 39 97-79

info@sukhamburg.com www.sukhamburg.com



# **LEGAL NOTICE**

## Copyright 2020 Schäfter+Kirchhoff GmbH. All rights reserved.

Text, image, graphic, sound, video and animation files and their arrangement on Schäfter+Kirchhoff GmbH webpages are protected by copyright and other protective laws. The content may not be copied for commercial use or reproduced, modified or used on other websites. [more]